


USB Free Link


DIRECT BALANCE – PC COMMUNICATION
NO ADDITIONAL PROGRAMS REQUIRED

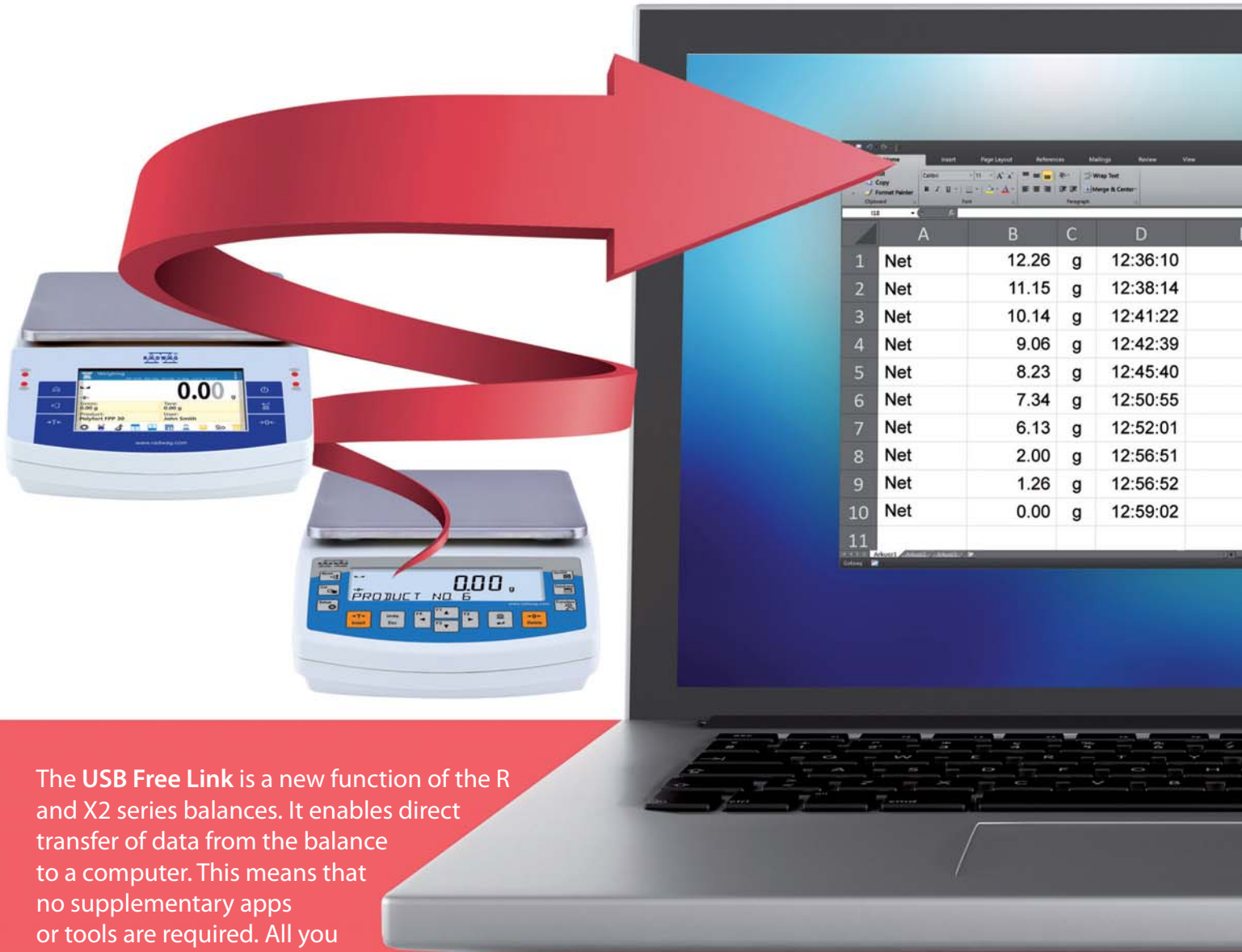

The **USB Free Link** is a new function of the R and X2 series balances. It enables direct transfer of data from the balance to a computer. This means that no supplementary apps or tools are required. All you need to do is simply to connect your balance and the computer using USB cable.

With the **USB Free Link** the data is transferred automatically to a currently operated computer app (e.g. Word, Excel, Notepad, etc.) upon pressing any balance key (e.g. Enter - printout, F1 - header, F3 - footer). The transferred text format remains unchanged.


Transfer of data to a currently operated application

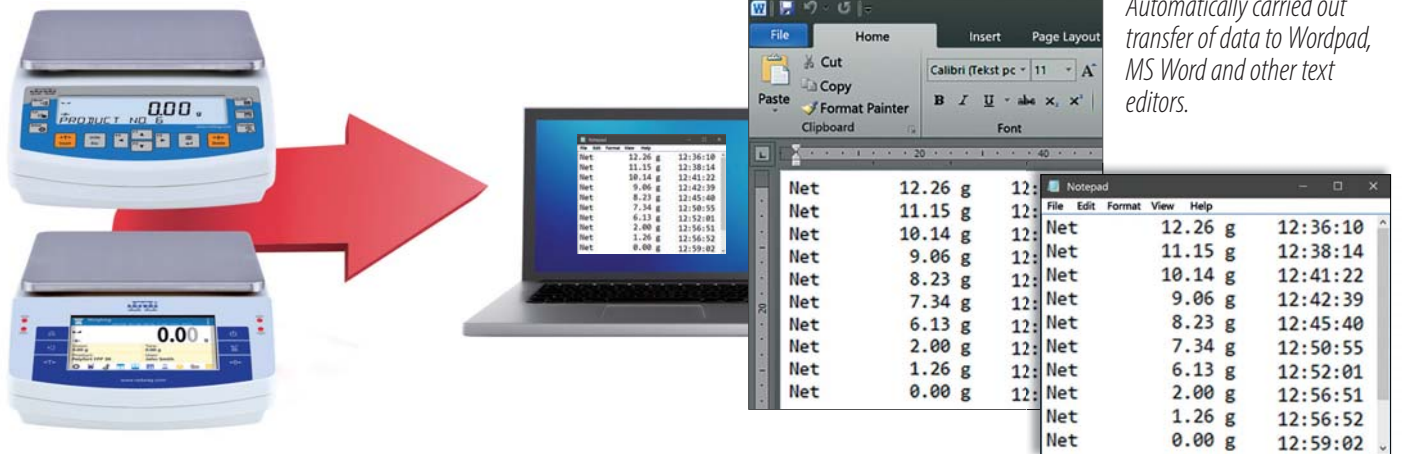

Automatically carried out transfer of data to Wordpad, MS Word and other text editors.

Net	12.26 g	12:36:10
Net	11.15 g	12:38:14
Net	10.14 g	12:41:22
Net	9.06 g	12:42:39
Net	8.23 g	12:45:40
Net	7.34 g	12:50:55
Net	6.13 g	12:52:01
Net	2.00 g	12:56:51
Net	1.26 g	12:56:52
Net	0.00 g	12:59:02

Transfer of data to a spreadsheet program

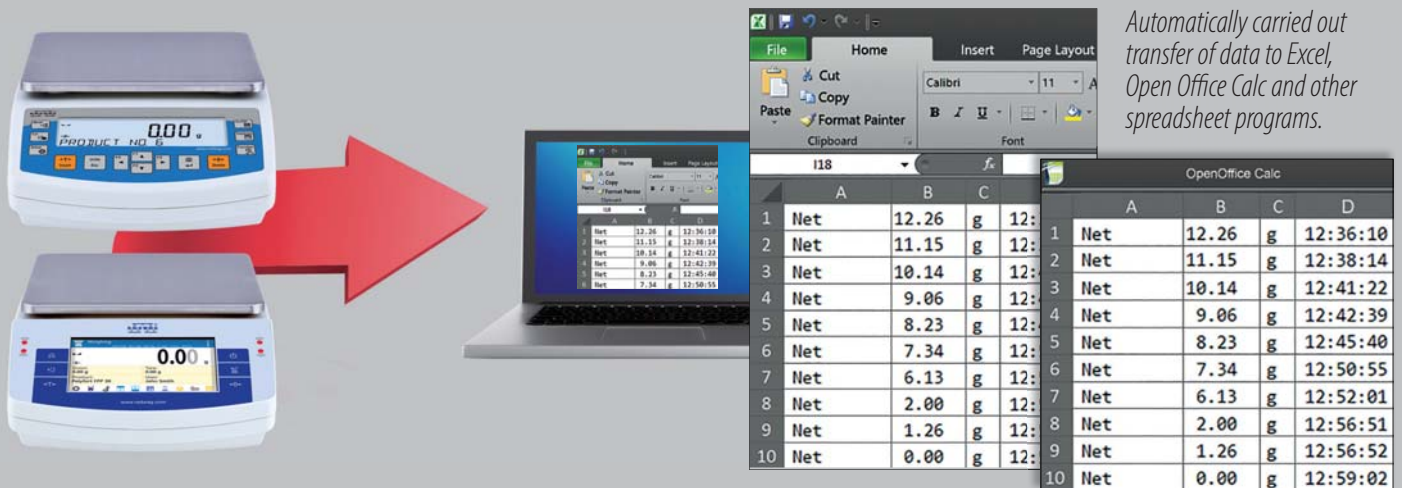

Automatically carried out transfer of data to Excel, Open Office Calc and other spreadsheet programs.

	A	B	C	D
1	Net	12.26 g	12:36:10	
2	Net	11.15 g	12:38:14	
3	Net	10.14 g	12:41:22	
4	Net	9.06 g	12:42:39	
5	Net	8.23 g	12:45:40	
6	Net	7.34 g	12:50:55	
7	Net	6.13 g	12:52:01	
8	Net	2.00 g	12:56:51	
9	Net	1.26 g	12:56:52	
10	Net	0.00 g	12:59:02	

Transfer of data to CSV and TXT files



Import of data to a freely selected spreadsheet program.